


TAYA 100 Series

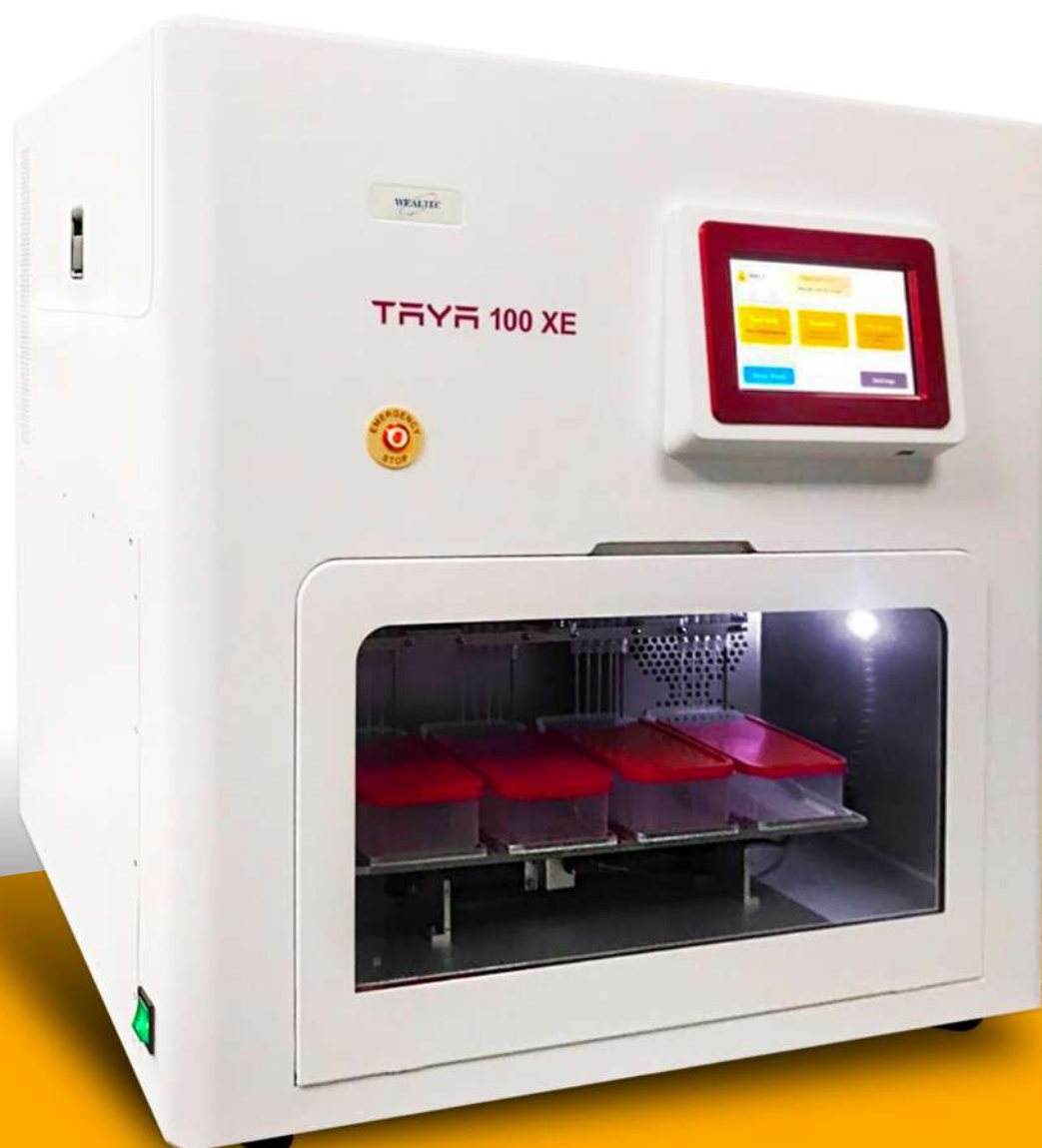

Automated Western Blot Processor

Automated Western Blot Processor

The TAYA 100 Series is an automated Western blot processor for incubating and washing membrane blots. The automation system simulates the conventional Western blotting procedures and allows users to customize their protocol settings, such as blocking and antibody incubation duration, washing duration and repeats, and options for recycling antibodies. If operator still prefer the conventional Western blotting procedures in their laboratories, the TAYA 100 Series preserves the techniques and provides consistent results with automation.



Increase Your Western Blot Quality

The TAYA 100 series simply brings your western blots to an impeccable quality. Bridging the gap between manual handling of dispensing and removing reagents from the membrane containers, the TAYA 100 series features fully programmable parameters for western blotting procedure.

Liberate Yourself from the Lab Bench

We are all aware of western blot is a tedious wet work. The TAYA 100 series automates reagents dispensing and removal. From the nitty-gritty of pouring and dumping reagents to monitoring the incubation time of sample blots, the TAYA 100 series ensures precise timing and liquid handling. The blots are ready for detection when you return to your lab bench.

Optional Cooling Chambers Model

The TAYA 100XE model features cooling chambers with temperature control for storing reagents and incubating membranes at 4°C for a prolong time. With this model letting a machine processing your Westerns is no longer a dream.









Features

Individual Probes

The individual probes effectively prevent cross-contamination and precisely dispensing, extracting, and recycling reagents, thanks to the diaphragm pumps and valve sets.

Blot container tray

The blot container tray is a programmable Two-Way Horizontal Shaking Device that can adjust the shaking speed and duration at each incubation or washing stage.

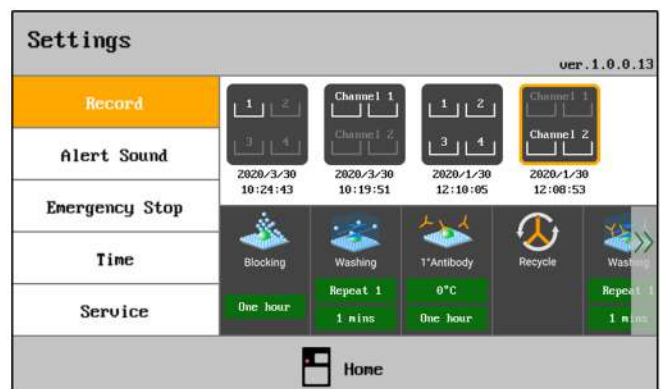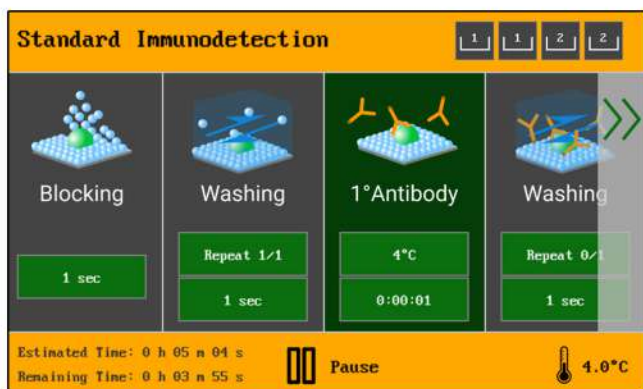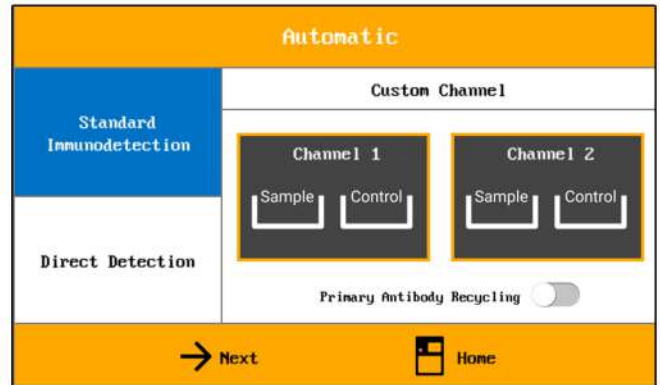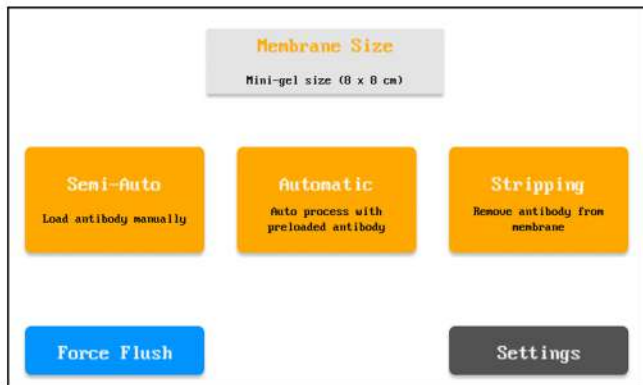

Easy-to-use User Interface

The TAYA 100 series comprises a built-in 7-inch LCD touchscreen computer, which features intuitive user interface and graphical western blotting layout. From the touchscreen, users can program the number of the membranes to use, shaking speed, number of washes, washing duration, and incubation time for each stage.



Specifications

| | TAYA 100XE | TAYA 100 |
|----------------------------------|-----------------------|------------------|
| Peltier controlled Temp. Module | 3 sets | N/A, R-T |
| Workflow mode selection | 10 workflow | 8 workflow |
| Control with touch panel | 7" LCD display | 7" LCD display |
| System condition display | Yes | Yes |
| Membrane pan | four | four |
| Primary Anti-Body add | Auto | Manual |
| Secondary Anti-body add | Auto | Manual |
| Primary Anti-body recollection | Yes | Yes |
| Stripping capability | Yes, 37°C | N/A |
| Auto waste drain | Yes | Yes |
| Adjustable shaking speed | Yes | Yes |
| Wash time adjustable | Yes | Yes |
| Cycles of wash adjustable | Yes | Yes |
| Temp. of reagent storage chamber | @4°C | RT |
| Reaction zone | 25°C / 0°C selectable | RT |
| Method save / recall | Yes | Yes |
| Emergency stop button | Yes | Yes |
| Dispensing channel | four | four |
| Waste drain channel | four | four |
| Operation Temperature | 0-45°C | 0-45°C |
| Physical dimensions (W*D*H) mm | 720 x 55 x 740 | |
| Power requirement | 96 - 240V, 200W | 96 - 240V / 120W |
| Weight | 65kg | 45kg |

The spec is subject to change without prior notice.

Ordering Information

| Catalog No. | Description |
|-------------|--|
| 1075001 | TAYA 100 XE, Automated Western Blot Processor, 96-240V, 50-60Hz. |
| 1075011 | TAYA 100 Western Blot Processor, 96-240V, 50-60Hz. |

Wealtec Corp.

1885 Meadowvale Way,
Sparks, NV 89431, USA

Tel: +1-775-351-2066 Fax: +1-775-351-2077

E-Mail:sales@wealtec.com

Wealtec Bioscience Co., Ltd.

27F., No.29-1, Sec. 2, Zhongzheng E. Rd.,
Danshui Dist., New Taipei City 251, Taiwan

Tel: +886-2-8809-8587 Fax: +886-2-8809-8589

Web-site:http://www.wealtec.com

